

**UNPACKING INSTRUCTIONS  
ILLUMINATOR BASE  
MODELS 7790-500 and 7790-502**

Upon receipt, carefully unpack the illuminator and retain all packaging material in the event the stirrer requires future shipment to Phipps & Bird for repairs.

Remove the protective paper sheet from the acrylic top.

The illuminator is shipped with the fluorescent light tube installed.

**INSTRUCTIONS FOR REPLACING FLUORESCENT LIGHT TUBE**

**DO NOT ATTEMPT TO INSTALL OR REMOVE LIGHT TUBE OR PERFORM ANY MAINTENANCE WHILE ILLUMINATOR IS CONNECTED TO A POWER SUPPLY.**

Turn illuminator upside down. Squeeze and lift white REFLECTOR SHIELD from the four TABS on illuminator FRAME. Turn LIGHT TUBE approximately one-quarter turn until end-pins of LIGHT TUBE align with slots on END BRACKETS. Slide the LIGHT TUBE from slots in END BRACKETS. Reverse the above procedure using a new replacement LIGHT TUBE. Replace reflector once LIGHT TUBE is installed.

**BE SURE TO READ THE NOTES ON PROPER ELECTRICAL GROUNDING.**

